IKATRON® ETC 1

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OPERATING INSTRUCTIONS GB 7

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D

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CE-DECLARATION OF CONFORMITY

GB

We declare under our sole responsibility that this product corresponds to the regulations 89/336EEC and 73/23EEC and conforms with the standards or standardized documents EN 61010-1; EN 50081; EN 50082; EN 55011 and EN 60555.

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Staufen, März 2002

Reiner Dietsche
Vice President

Wolfgang Buchmann
Corporate Director - Quality Assurance

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Guarantee

You have purchased an original IKA laboratory machine which meets the highest engineering and quality standards.

In accordance with IKA guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

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
Correct use

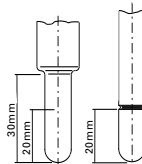The IKATRON ETC 1 is an instrument for the precise control of temperatures. The sensor of the instrument must be immersed at least 20mm deep in the medium to be tempered. The device can only be connected to special IKA magnetic stirrers. The magnetic stirrer must be approved for connecting the IKATRON ETC 1 (for example RH KT - RH KT/C)

Unpacking

Please unpack the machine carefully and inspect for damage. It is important that any transit damage should be noted at the time of unpacking. In certain circumstances it may be necessary to investigate immediately (post, rail or freight forwarder).

As supplied, the package contains: one IKATRON ETC 1 and one set of operating instructions.

Commissioning

Plug the instrument into the contact thermometer jack on the magnetic agitator (e.g. IKA RH KT - RH KT/C). The instrument is activated when the agitator is switched on.

An automatic self-test is performed when the instrument is activated. During the self-test all the segments of the LCD display are visible (**Fig.1**). Then the LCD displays illustrated in **Fig.2** to **Fig.3** appear (standard factory setting).

Useful information

The IKATRON ETC 1 works as a two-point controller with a small hysteresis. In order to achieve good results, please note the following points:

- Use vessels with flat bottoms.
- The heating temperature of the magnetic stirrer should be adjusted to about twice the setpoint of the IKATRON ETC 1.
- For rapid heating up, or if the setpoint temperature is not being reached, the temperature of the heating plate should be increased.
- To avoid exceeding the setpoint temperature initially, or to prevent the temperature from moving back and forth between too high and too low, the temperature of the heating plate should be reduced.
- When controlling the temperature of oils, vessels with optimal properties for conducting heat should be used, for example IKA's H15.
- If possible, the viscosity of the oils should be less than 50mPas, and they should be stirred at high speeds (to ensure adequate mixing).

Setting the nominal temperature

The nominal temperature is adjusted in 1K steps by pressing the "+" and "-" keys. If the key is held down for longer than 5 seconds, the nominal temperature setting changes in 10K steps. The LCD display shows: the

current actual temperature with physical units, the nominal temperature at top right, and OFF to indicate that the heater of the connected device is switched off. (Fig. 5)

If the set nominal temperature is higher than the current actual temperature, the word OFF in the display changes to ON and the heater of the connected device is switched on. (Fig. 4)

Replacing sensor

In order to replace the sensor, pull the sensor and the cap downward until the plastic snap fastener is released. Then push the sensor with the cap over the fitting on the instrument to reestablish the connection.

Error messages

The following error messages may occur (Fig. 6/7) and have the causes listed below:

- Er 1 Sensor not connected or defective (open circuit).
 - Er 3 Internal temperature above +76 °C.
 - Er 6 Sensor short circuit.
 - Er 8 Calibration error (re-calibration by field service required).
 - Er 9 IC107 (4066) defective
 - Er 11 Supply voltage <7.5V
 - Er 12 Supply voltage >16.1V
- not comp the instrument is not compatible

Accessories

- H 62 ETC 1** Standard sensor 270 mm stainless steel
- H 66 ETC 1** Glass sensor 270 mm glass

Technical data

Measuring range:	°C	0 ... 350
Resolution:	K	0.1 (0 ... +100 °C)
	K	1 (100 ... 350 °C)
Measuring accuracy:	K	±0.3 + sensor tolerance Pt 1000 DIN IEC 751 Class B
Setting accuracy:	K	1
Control deviation:	K	± 3
Supply voltage:	V DC	8 ... 16
Current consumption:	mA	16 (at 9V)
Ambient temperature:	°C	0 ... 60
Relative humidity:	%	80
ON time:	%	100
Degree of protection:		IP 44
Cable length:	mm	350
Plug:		6-pin DIN 45 322
Connection:		on special IKA magnetic stirrer
Dimensions: (W x D x H)	mm	70 x 65 x 150 (without sensor)
Weight:	kg	0.2

List of spare parts

Item	Designation	Item	Designation
1	Rear wall	13	Plastite screw 2.5 x 8
2	Clamp	14	H 62 ETC 1 stainless steel sensor Pt1000
3	Sealing ring	15	Cap
4	Spiral cable	16	Gasket
5	Sealing insert		
6	Union nut		
7	Housing		
8	Dished washer		
9	Membrane keypad		
10	Sensor jack; complete		
12	BLP temperature regulator		

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